

Product Information

Pepsin from porcine gastric mucosa

Lyophilized powder

P7012

Product Description

CAS Registry Number: 9001-75-6

Enzyme Commission (EC) Number: 3.4.23.1

Molecular Mass: 34,620 (calculated from amino acid sequence)¹pI:^{2,3} 2.2-3.0; 2.2, 2.8 λ_{max} : 278 nm⁴Extinction coefficient: $E^{\text{M}} = 51,300^4$

Synonym: Pepsin A

Pepsin, unlike some other peptidases, hydrolyzes only peptide bonds, not amide or ester linkages. The cleavage specificity includes peptides with an aromatic acid on either side of the peptide bond, especially if the other residue is also aromatic or a dicarboxylic amino acid. Increased susceptibility to hydrolysis occurs if there is a sulfur-containing amino acid close to the peptide bond, which has an aromatic amino acid. Pepsin will also preferentially cleave at the carboxyl side of Phe and Leu, and to a lesser extent at the carboxyl side of Glu residues. Pepsin will not cleave at Val, Ala, or Gly linkages.⁵ Some good substrates of pepsin include:

- Z-L-tyrosyl-L-phenylalanine
- Z-L-glutamyl-L-tyrosine
- Z-L-methionyl-L-tyrosine

Amidation of the C-terminal carboxyl group prevents hydrolysis by pepsin.^{5,6}

Pepsin is commonly used in the preparation of Fab fragments from antibodies. The optimal pH for the pepsin reaction is 1.5-2.5, which will not be detrimental to the antibody if it is not exposed for long durations to the low pH. Solutions should be adjusted to neutral pH for storage. The control of pepsin digestion of antibodies has been reported.⁷

For general protein digestion, suggested conditions are a 0.4% solution of pepsin in 10 mM HCl, and digestion for 30-90 minutes at 37 °C. Pepsin has an optimal activity with native proteins at pH ~1.0. However, with some denatured proteins, the optimal activity is in the range of pH 1.5-3.5.^{8,9} Pepsin is inhibited by several Phe-containing peptides.¹⁰

Several theses¹¹⁻¹⁵ and dissertations¹⁶⁻²² have cited use of product P7012 in their protocols.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

Pepsin is soluble in deionized water at 10 mg/mL, and 4 mg/mL in cold 10 mM HCl.

Storage/Stability

In lyophilized form as sold, this product should be stored at -20 °C.

Solutions of pepsin at pH 4.4 are stable at -20 °C for about 2-3 months.²³ Pepsin is not active when not at an acidic pH and a solution is stable at pH 6-7. Bringing the pH up to 8, however, will irreversibly inactivate pepsin. Pepsin is irreversibly denatured at pH 8.5-11 at room temperature.²⁴

References

1. Sepulveda, P. *et al.*, *J. Biol. Chem.*, **250**(13), 5082-5088 (1975).
2. Jonsson, M., *Acta Chem. Scand.*, **26**(9), 3435-3440 (1972).
3. Malamud, D., and Drysdale, J.W., *Anal. Biochem.*, **86**(2), 620-647 (1978).

4. Perlmann, G.E., *Proc. Natl. Acad. Sci. USA*, **45(7)**, 915-922 (1959).
5. Sweeney, P.J., and Walker, J.M., *Methods Mol. Biol.*, **16**, 271-276 (1993).
6. Dixon, M., and Webb, E.C. (eds.), *Enzymes*. Academic Press (New York, NY), p. 262 (1979).
7. Rea, D.W., and Ultee, M.E., *J. Immunol. Methods*, **157(1-2)**, 165-173 (1993).
8. Christensen, L.K., *Arch. Biochem. Biophys.*, **57(1)**, 163-173 (1955).
9. Schlamowitz, M., and Peterson, L.U., *J. Biol. Chem.*, **234(12)**, 3137-3145 (1959).
10. Knowles, J.R. *et al.*, *Biochem. J.*, **113(2)**, 343-351 (1969).
11. Gangloff, Mary Beth, "Composition and *in vitro* Digestion of Barley, Oat, and Wheat Brans". Oregon State University, M.S. thesis, p. 69 (1992).
12. Karava, Nilesh B., "The effect of heating chicken muscle on formation of bioavailable forms of iron". University of Massachusetts Amherst, M.S. thesis, pp. 29, 47, 70, 94 (2008).
13. Robinson, Mary Anna, "Production, Fractionation, and Evaluation of Antioxidant Potential of Peptides Derived from Soy Protein Digests". University of Waterloo, M.App.Sc. thesis, pp. 22, 84, 85, 111 (2010).
14. Kodali, Mahesh Chandra, "Identification and Quantification of the Collagen Type I, III and V in Rabbit Patellar Tendons". Wright State University, M.S. thesis, p. 34 (2015).
15. Bhoi, Rahulkumar Kantibhai, "Characterization & extraction of extracellular matrix from porcine adipose tissue". University of Texas at El Paso, M.Eng. thesis, p. 16 (2017).
16. Ahn, Joomi, "Local hydrogen deuterium exchange mass spectrometry: from pressurized online digestion to pepsin proteolysis". Northeastern University, Ph.D. dissertation, p. 157 (2012).
17. Helbig, Anne, "Digestion of dietary fat: Gastrointestinal behaviour of emulsions and human physiological responses". Wageningen University, Ph.D. dissertation, p. 51 (2013).
18. Marcolini, Elena, "NMR based foodomics to investigate the digestibility of protein-rich food products". Università di Bologna, Ph.D. dissertation, p. 68 (2015).
19. Alhasawi, Fatemah M., "Influence of food matrix physico-chemical properties on the biophysics of digestion and *in situ* luminal viscosity". Rutgers, the State University of New Jersey, Ph.D. dissertation, pp. 67, 71, 95, 120 (2018).
20. Molina, Catalina Pineda, "Characterization of the host tissue response induced by a biosynthetic material composed of poly (4-hydroxybutyrate)". University of Pittsburgh, Ph.D. dissertation, p. 18 (2018).
21. Lee, Jung Wook, "Optimization of Canola Co-Product Utilization in Swine". South Dakota State University, Ph.D. dissertation, p. 48 (2019).
22. Kellaway, Simon Christopher, "Hydrogels derived from decellularised tissues for nerve repair". University College London, Ph.D. dissertation, p. 72 (2021).
23. Rajagopalan, T.G. *et al.*, *J. Biol. Chem.*, **241(21)**, 4940-4950 (1966).
24. Ryle, A.P., *Methods Enzymol.*, **19**, 316-336 (1970).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.
P7012pis Rev 04/22 ARO,RXR,GCY

MERCK